

Aug 14, 2024

Version 1

Wheat spike meristem micro-dissection V.1

DOI

<https://dx.doi.org/10.17504/protocols.io.3byl49r2zgo5/v1>



Isabel Faci¹, Anna Elisabeth Backhaus², Cristobal Uauy¹

¹John Innes Centre, Norwich Research Park, Norwich NR4 7UH, United Kingdom; ²ICARDA

Uauy Lab - JIC



Isabel Faci

John Innes Centre

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.3byl49r2zgo5/v1>

Protocol Citation: Isabel Faci, Anna Elisabeth Backhaus, Cristobal Uauy 2024. Wheat spike meristem micro-dissection. protocols.io <https://dx.doi.org/10.17504/protocols.io.3byl49r2zgo5/v1>



License: This is an open access protocol distributed under the terms of the **[Creative Commons Attribution License](#)**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: July 30, 2024

Last Modified: August 14, 2024

Protocol Integer ID: 104289

Keywords: Meristems, Early development, Wheat, Dissection, developing meristem, wheat, spike, barley, growing leaf, spikes during early stage

Funders Acknowledgements:

UK Biotechnology and Biological Sciences Research Council (BBSRC)

Grant ID: BB/X011003/1

UK Biotechnology and Biological Sciences Research Council (BBSRC)

Grant ID: BB/X01102X/1

European Research Council

Grant ID: 866328

Disclaimer

This is one of many ways to dissect wheat (inflorescence) meristems. Please note that genotypes might differ in appearance from the figures provided. We hope this protocol provides a starting point which can then be adapted as needed by individual labs.

Abstract

The aim of this protocol is to micro-dissect young wheat plants to access developing meristems which are found enveloped within growing leaves. This protocol can be used to dissect spikes during early stages of development, for both wheat and barley.



Materials

- Dissection knives: GeeEdge ophthalmic slit knife model CJY-01-2.2. Another recommended knife is the GeeEdge MVR CJY-04-20g(Z), up to personal preference. We re-use them until they break. They can be purchased in AliExpress.
- Stereo microscope (with additional lighting and camera): We use Leica S9 (with 0.6X, 1X, 2X, 3X, 4X, 5.5X) with a HXCAM HiChrome HR4 Lite camera and a Photonic Optics light source.
- Software GX-Capture, to take pictures.
- Small pair of scissors.
- Optional (if doing #step 7): RNase removal BLITZ
- Optional (if doing #step 7): metallic beads, previously cleaned with 0.5 M HCl and 70% ethanol. Two per rep, either 3 mm or 5 mm diameter (*Hardened 52100 Chrome Steel Ball Bearings SIMPLY BEARINGS LTD*)
- Optional (if doing #step 7): liquid nitrogen
- Optional (if doing #step 7): dry ice

Abstract

- 1 Cereal plant development and growth rate are affected by environment, genetics and their interaction. A plethora of decisions are made in early stages of inflorescence/spike development, when the meristem primordium is being differentiated into spikelet and inflorescence meristems, depending on the environmental cues and genetics (Kirby, 1984). To study the factors that affect these developmental decisions, we usually micro-dissect, in-depth phenotype and/or collect spike meristems at very early developmental stages (*Note 1*). This protocol (adapted from Kirby, 1984) can be used to dissect spikes during early stages of development (up to approximately Waddington scale stage 4.0., (Waddington et al., 1983))(*Note 2*). It will vary slightly vary between stages and we will specify these differences throughout the protocol.

Note 1: The dissected spike meristems can be collected and pooled into cool (dry-ice) 2 mL Eppendorf tubes to perform RNA extraction or nuclei isolation (briefly described in Step #7). This specific protocol does not describe the RNA extraction or nuclei isolation.

Note 2: Due to the high similarity in wheat and barley inflorescences, this protocol would also be suitable for barley.

Waddington stages:  Waddington-scale.tif 1.7MB

Cereal developmental guide (Kirby, 1984):

 Cereal Development Guide KIRBY 1... 15.8MB

Materials needed - Setup

- 2 Gather the following materials:
 - **Dissection knives:** GeeEdge ophtalmic slit knife model CJY-01-2.2. Another recommended knife is the GeeEdge MVR CJY-04-20g(Z), up to personal preference. *These are one-use knives, however we clean and re-use them until the blade is worn off (they last for months). They can be purchased in AliExpress.*
 - **Stereo microscope** (with additional lighting and camera): We use Leica S9 (0.6X, 1X, 2X, 3X, 4X, 5.5X) with a HXCAM HiChrome HR4 Lite camera and a Photonic Optics light source.
 - **Software GX-Capture:** to take pictures.
 - Small pair of **scissors**.
 - **RNase removal** spray (BLITZ) (*Optional*): Only needed if meristems will be collected for RNA extraction, nuclei isolation, or similar (explained in #7). If used, make sure to

use it to clean all your surfaces and knives.

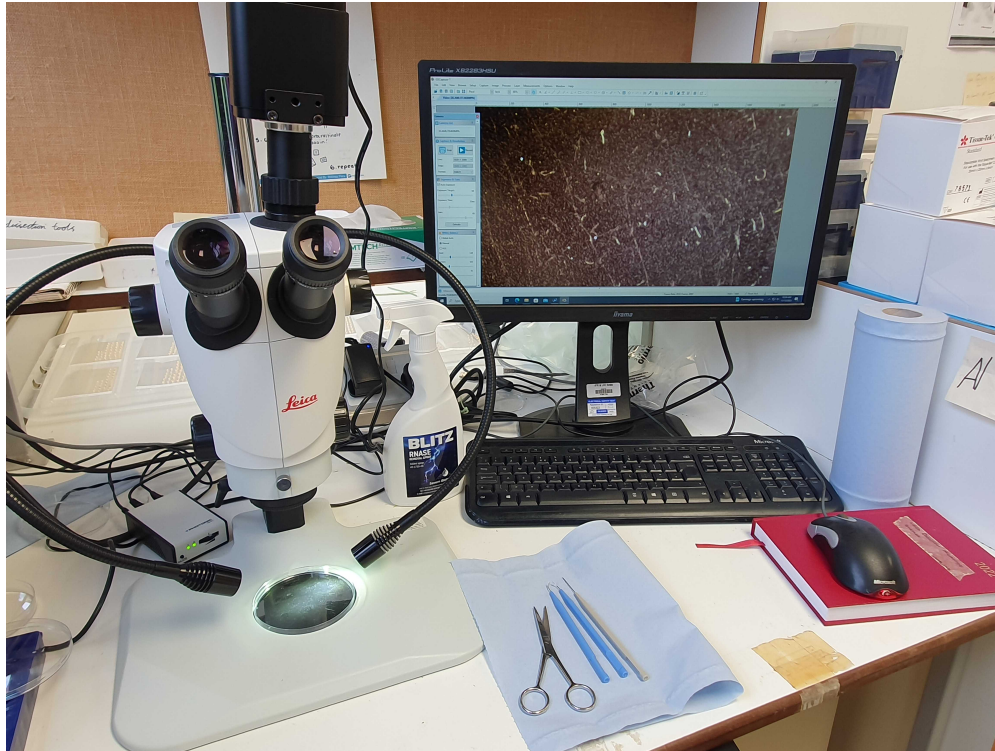


Fig. 1. Set up: stereo microscope, scissors, dissection knives, PC with GX-Capture software.

Dissection

- 3 Examine plant and **identify the main tiller. Cut just below roots** (Fig. 2). It is important to leave some roots to be able to hold the plant in later steps under the microscope. Clean the roots with some tissue to avoid having soil on the work area.



Fig. 2. Single main tiller with cut roots.

- 4 **Remove leaves one by one, gently.** The aim is to peel back and remove leaves so you can access the developing meristem which is within them. You can do this with the knife (CJY-01-2.2), and helping with your hands. Always make sure that the highest node (the node furthest away from the roots) is still there and intact, as this is where the spike meristem is (*Note 3*). In early stages of development (approx. up to W2.5-3), the highest node is still at the very base of the tiller. In later stages, this node will start to go up (as observed in Fig. 3).

Note 3: You can often feel with your fingers where the nodes are, but this is not always the case. For example, in Fig. 3, it is easy to feel the node indicated with orange, but not the one indicated in white. Because of this, it is recommended to practise this technique and to always be careful with the whole tiller until you find the highest node.

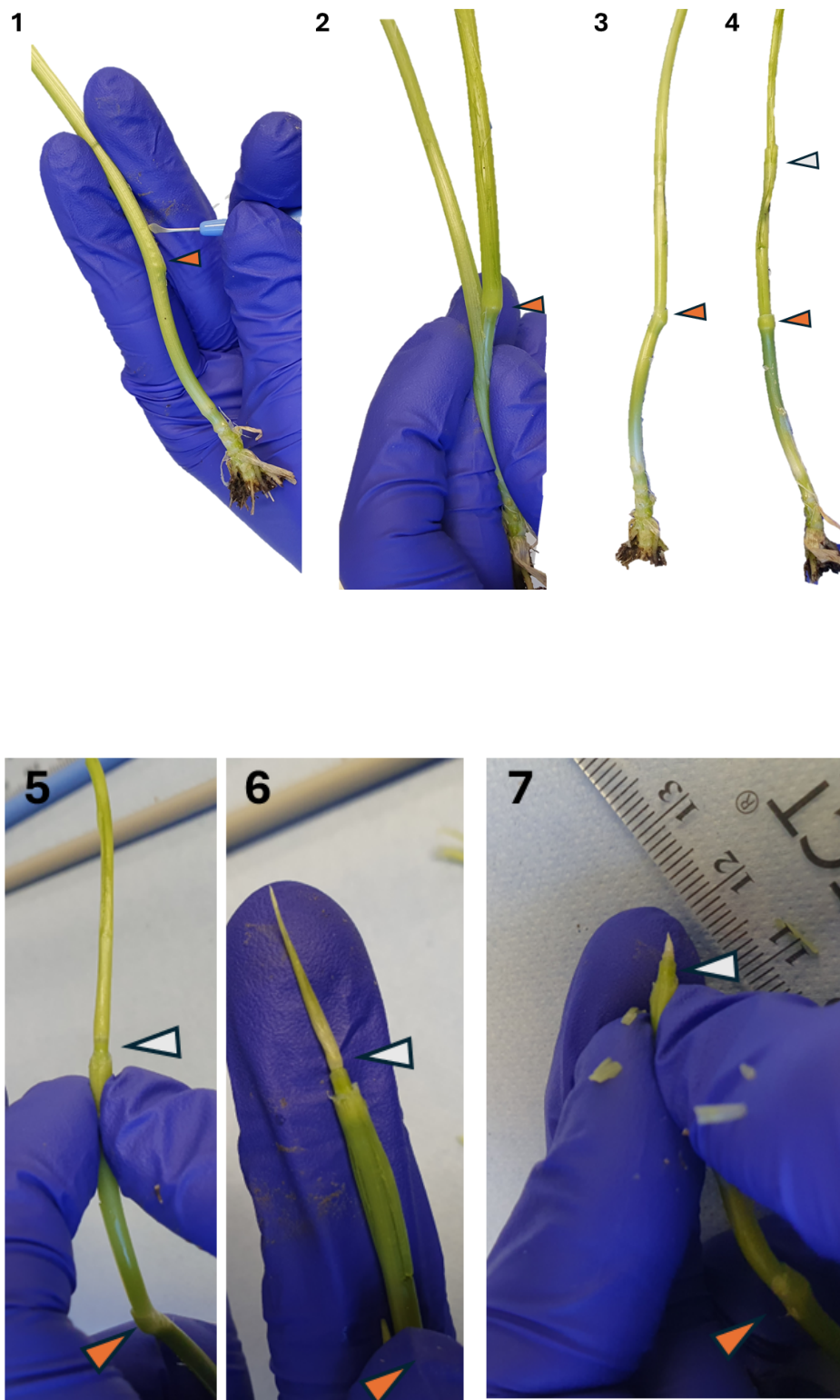


Fig. 3. Removing leaves to access meristem. Triangles mark nodes. The spike meristem is just above the white triangle. In step 7, there are still a couple of intact leaves which envelop the meristem and which will be removed carefully under the microscope.

Once you remove most leaves and the next leaf is ~ 2 mm (step 7 in Fig. 3), continue dissecting under the stereo microscope.

- 5 **Dissect under the microscope.** Use a Petri dish to place the sample and hold the roots/stem with one hand. Hold the knife with the other hand (Fig. 4), and dissect carefully, like in the videos below.

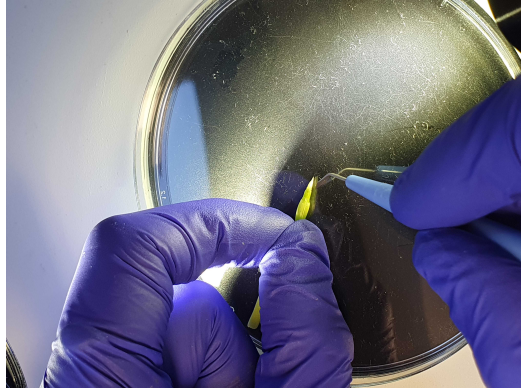


Fig. 4. Dissection under the microscope.

To see an example of how to dissect under the microscope, **watch these videos**, taken with the microscope camera:

https://www.youtube.com/embed/vZ784Dj9cZY?si=7tB786BuwGlg5-_q

<https://www.youtube.com/embed/XQ06DrLibro?si=b3Zj9ITCdvZQ-Cyp>

- 6 **Capture images with the camera.** These are examples of pictures of the developing meristems:

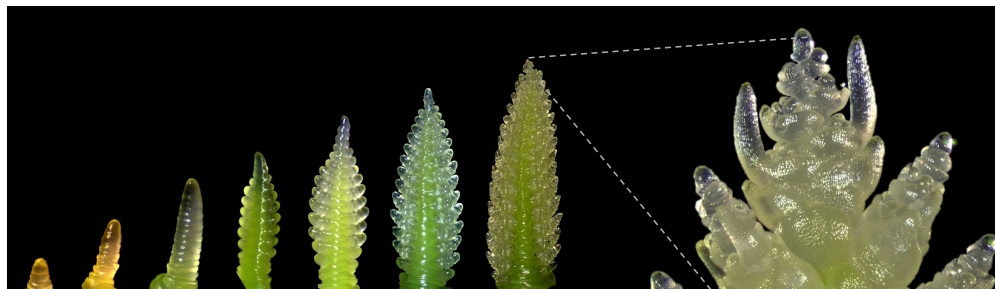


Fig. 5. Dissected meristems captured with GX-Capture under Leica S9 stereo microscope.

Collection in clean and cold conditions (optional)

- 7 To collect for RNA extraction or similar, make sure that the knives used and all surfaces in the setup are previously cleaned with **RNase removal**. Clean the surfaces and knives between the collection of each meristem.



Once the meristem is visible, like in the video, cut the spike at the base with a clean knife. Then, hold a RNase-free small pipette tip and pick up the meristem to place it carefully into cool (dry-ice) labelled 2 mL RNase-free Eppendorf tubes to perform RNA extraction, nuclei isolation or equivalent molecular biology processing of the sample.

For these protocols, we often pool several meristems at the same stage into a tube. We also place two nitrogen-cooled metallic beads (previously cleaned with 0.5 M HCl and 70% ethanol) in each tube. The reason why we add two beads to each tube is because the first step of RNA extraction or nuclei isolation will be grinding the tissue (in cold conditions). The metallic beads are the method that works best for us for grinding low amounts of wheat meristems.

Once a sample is collected, we place it in liquid nitrogen. The sample can be processed or long-term stored in -80 °C, until processed.

This specific protocol does not describe the RNA extraction or nuclei isolation.

Protocol references

Kirby, E.J.M., 1984. *Cereal development guide*, 2nd edition. Stoneleigh, Kenilworth, UK, NAC Cereal Unit. 85pp.