


Jun 24, 2025

🌐 Whole Genome Sequencing of Hepatitis B Virus

DOI

<https://dx.doi.org/10.17504/protocols.io.e6nvwx1zvmk/v1>

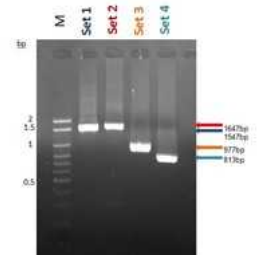

Myat Htut Nyunt¹, Wah Wah Aung², Yi Yi Kyaw²

¹Advanced Molecular Research Centre, Department of Medical Research, Republic of the Union of Myanmar;

²Department of Medical Research



MYAT NYUNT

DMR

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.e6nvwx1zvmk/v1>

Protocol Citation: Myat Htut Nyunt, Wah Wah Aung, Yi Yi Kyaw 2025. Whole Genome Sequencing of Hepatitis B Virus. protocols.io <https://dx.doi.org/10.17504/protocols.io.e6nvwx1zvmk/v1>

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited



Protocol status: Working

We use this protocol and it's working

Created: June 24, 2025

Last Modified: June 24, 2025

Protocol Integer ID: 220786

Keywords: HBV, whole genome sequencing, hepatitis B virus, next-generation sequencing, sequencing, hbv across different genotype, universal primers for hbv, procedure for whole genome sequencing, standardized method for whole genome sequencing, generation sequencer, hbv, genome, different genotype, genotype, using universal primer

Abstract

This standard operating procedure (SOP) describes the procedure for whole genome sequencing (WGS) of the hepatitis B virus using a next-generation sequencer. Next-generation sequencing (NGS) has been widely used for genome-wide studies in many aspects of infectious as well as non-communicable diseases. However, there is no standardized method for whole genome sequencing of HBV across different genotypes. This SOP was developed using universal primers for HBV genomes across genotypes A to G.

Guidelines

Follow the Good Laboratory Practice in all steps.

Use powder-free gloves to handle the reagents and all libraries.

Use filtered-tips for pipetting.



Materials

1. General

- Micropipettes and tips (10 μ L, 200 μ L, 1000 μ L)
- Disposable powder-free gloves
- 2N NaOH stock
- Microcentrifuge tubes (low bind)
- 15ml tube
- Vortex
- Microcentrifuge
- Centrifuge
- PCR plates with adhesive seals
- Kimwipes
- Lens tissue

2. Targeted amplifications

- GoTaq Long PCR Master Mix (M4021)
- Primers
 - Set 1: 251F (GAC TYG TGG TGG ACT TCT C)
1797R (CCA ATT TMT GCY TAC AGC CTC)
 - Set 2: 2300F (CCA CMW AAT GCC CCT ATC)
654R (GSC CCA MBC CCA TAG G)
 - Set 3: 1859F (ACT NTT CAA GCC TCC RAG CTG)
2835R (GTT CCC AVG WAT AWG GTG AYC C)
 - Set 4: 1584F (ACT TCG MBT CAC CTC TGC ACG T)
2396R (GTC KGC GAG GYG AGG GAG TT)

3. Library Preparation Kit

- Illumina Microbial Amplicon Prep (iMAP)
- PhiX Control Kit, version 3 (10 μ L at 10 nM) (#15017666)
- Illumina DNA/RNA UD Indexes Sets A–D (Catalog nos. 20091654, 20091656, 20091658, and 20091660)
- MiSeq Reagent Kit, version 2 (300 cycles)

4. Other reagents

- Squeeze bottles
- pH sticks
- Molecular grade distilled water
- Tween 20
- Isopropyl Alcohol
- Absolute Ethanol
- Ice bucket with ice
- Water bath

Before start

DNA extraction should be done accordingly using the QIAamp DNA Kit (QIAGEN) in accordance with the manufacturer's instructions.



General Precaution before the procedure

- 1 Follow the Good Laboratory Practice in all steps.
- 2 Use powder-free gloves to handle the reagents and all libraries.
- 3 Use filtered-pipette tips for pipetting.
- 4 DNA extraction should be done accordingly using the QIAamp DNA Kit (QIAGEN) in accordance with the manufacturer's instructions.

Amplification

- 5 PCR preparation
 - 5.1 Thaw the Long PCR mastermix.
 - 5.2 Prepare the PCR reactions as follows: Long PCR mix (2X), DNA template, DDW, Total 20 μ L.
- 6 PCR Amplification
 - 6.1 Prepare the PCR conditions as follows: Initial denaturing 02:00, 01:00, 01:00 35 cycles, 01:15, 05:00, Hold.
- 7 Gel-electrophoresis
 - 7.1 Confirm the amplification of the samples using 1% agarose gel-electrophoresis.

Library preparation

8 Tagment PCR amplicons

- 8.1 5 μ L each from four PCR amplicons were pooled together in a tube. (Total 20 μ L)
- 8.2 Thaw EBLTS HT (4°C) at RT and vortex thoroughly before use.
- 8.3 Thaw TB1HT (-20°C) at RT and vortex thoroughly before use.
- 8.4 Prepare the reagent mix: 12 μ L TB1 HT, 4 μ L EBLTS HT, 20 μ L H₂O.
- 8.5 Use 30 μ L of reagent mix to each sample (Add 30 μ L of mix to the sample) (Total 50 μ L).
- 8.6 Seal and shake 1600 rpm for 1 min.
- 8.7 Centrifuge 1000 g for 1 min.
- 8.8 Run: 55°C for 5 min, 10°C Hold (lid-heated, 50 μ L volume).

9 Post Tagmentation Cleanup

- 9.1 Keep ST2 HT (RT) vortex before use.
- 9.2 TWB HT (4°C) at RT and vortex before use.
- 9.3 Centrifuge the plate for 1 min.
- 9.4 Add 10 μ L of ST2 HT.

- 9.5 Seal and shake.
- 9.6 Centrifuge.
- 9.7 Incubate at RT for 5 min.
- 9.8 Place on magnetic stand until clear (~3 min).
- 9.9 Remove supernatant.
- 9.10 Wash the bead by: Remove from magnetic stand, Add 100 μ L TWB, Seal and shake (pipette mix), Centrifuge prn, Place on magnetic stand until clear (about 3 min), Remove supernatant.
- 9.11 Wash again a second time, but leave the supernatant until the next step.
- 10 Amplify Tagmented Amplicons
- 10.1 Thaw EPM HT (-20°C) on ice until use (mix by inverting).
- 10.2 Index 10 bp adapter (-20°C), thaw RT and centrifuge after vortex.
- 10.3 Prepare the PCR mix by 20 μ L of EPM + 20 μ L of H₂O.
- 10.4 Vortex to mix the PCR mix.
- 10.5 Remove the supernatant wash in previous cleanup.

- 10.6 Use 10 μ L tips to remove residual wash solution.
- 10.7 Remove from magnetic stand.
- 10.8 Add 40 μ L of PCR mix.
- 10.9 Add 10 μ L of index.
- 10.10 Seal and shake.
- 10.11 Centrifuge briefly.
- 10.12 Run inside a thermocycler as follows: 72°C for 3 min, 98°C for 3 min, 7 cycles of 98°C for 20 sec, 60°C for 30 sec, 72°C for 1 min, 72°C for 3 min, Hold at 10°C (lid heated 100°C, vol 50 μ L).
- 11 Pool and cleanup libraries
 - 11.1 Take 7 μ L each from libraries and pool to 1.5 ml tube.
 - 11.2 Add 0.9 x of the IPB.
 - 11.3 Shake well and briefly spin down.
 - 11.4 Incubate at RT for 5 min.
 - 11.5 Place on magnetic stand for 3 min.



- 11.6 Discard supernatant.
- 11.7 Add 500 μ L of freshly prepared 80% Ethanol (while keeping the tube on magnetic stand).
- 11.8 Remove the supernatants.
- 11.9 Add a second time 500 μ L of 80% Ethanol.
- 11.10 Remove the supernatant.
- 11.11 Use 10 μ L pipette to remove extra reagents.
- 11.12 Air dry (about 3 min).
- 11.13 Add RSB 55 μ L to the bead after removing from magnetic stand.
- 11.14 Incubate 2 min.
- 11.15 Place to magnetic stand for 2 min.
- 11.16 Elute 50 μ L of pooled libraries.
- 11.17 Check Qubit and the initial result showed too high. So 5 μ L of the library was diluted with 25 μ L of the EBT and measured again.
- 11.18 After diluting to 4 nM, recheck the concentration. (ng/ μ L that is equal to nM. We proceed accordingly).
- 12 Denaturing and running in MiSeq



- 12.1 Prepare the 0.2N NaOH (0.2N NaOH = 100 μ L of 1N NaOH + 400 μ L of LGW).
- 12.2 5 μ L of lib + 5 μ L of 0.2N NaOH.
- 12.3 Incubate for 5 min at RT.
- 12.4 Add prechilled HT 990 μ L (to become 24 pM, 1000 μ L).
- 12.5 Combine 311 μ L of 24 pM + 288 μ L of prechilled HT (600 μ L of 12.5 pM lib).
- 12.6 Prepare 12.5 pm (100 μ L) of PhiX (by dilute 63 μ L of 20 pM + 38 μ L of HT).
- 12.7 Add 5% (30 μ L) of PhiX control to load.

Protocol references

MiSeq System: Denature and Dilute Libraries Guide. Illumina Document #15039740 v01.

Chook JB, Teo WL, Ngeow YF, Tee KK, Ng KP, Mohamed R. Universal primers for detection and sequencing of hepatitis B virus genomes across genotypes A to G. J Clin Microbiol. 2015 Jun;53(6):1831–1835. doi:10.1128/JCM.03449-14

Illumina, Inc. Protocol (v00): Illumina Microbial Amplicon Prep (IMAP) library preparation protocol [Internet]. San Diego (CA): Illumina; 2023 [cited 2025 Jun 16]. Document No. 200039808.
Available from: https://support-docs.illumina.com/LP/IMAP/Content/LP/IMAP/Protocol%28LibPrep%29_IMAP.htm