

Jan 07, 2019

Version 1

Wholemount Immunolabeling for GUT Samples V.1

DOI

<https://dx.doi.org/10.17504/protocols.io.wr6fd9e>

Marthe Howard¹

¹University of Toledo Health Sciences Campus



Marthe Howard

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.wr6fd9e>

Protocol Citation: Marthe Howard 2019. Wholemount Immunolabeling for GUT Samples. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.wr6fd9e>

Manuscript citation:

Development. 2011 Nov;138(21):4789-800. doi: 10.1242/dev.060053. Targeted deletion of Hand2 in enteric neural precursor cells affects its functions in neurogenesis, neurotransmitter specification and gangliogenesis, causing functional aganglionosis. Lei J1, Howard MJ.

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: January 04, 2019

Last Modified: January 07, 2019

Protocol Integer ID: 18974

Keywords: wholemount immunolabeling, enteric explants, wholemount immunolabeling for gut sample, immunostaining large block, mouse gut tissue, large pieces of mouse gut tissue, gut sample, large blocks of tissue, gut, tissue, tissues shold

Abstract

Wholemount Immunolabeling-Gut

This protocol is used for immunostaining large blocks of tissue. The application that it was developed for is large pieces of mouse gut tissue. Tissues shold be well washed prior to fixation.

Attachments



Wholemount
immunosti...

220KB

