



Jan 09, 2023

Wisecaver Lab DNAzol-based extraction of high molecular weight DNA from photosynthetic sea slugs

DOI

dx.doi.org/10.17504/protocols.io.n92ldprnl5b/v1

Jennifer H Wisecaver¹, Raeya Ogas¹

¹Purdue University

Wisecaver Lab



Jennifer H Wisecaver

Purdue University

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.n92ldprnl5b/v1>

Protocol Citation: Jennifer H Wisecaver, Raeya Ogas 2023. Wisecaver Lab DNAzol-based extraction of high molecular weight DNA from photosynthetic sea slugs. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.n92ldprnl5b/v1>

License: This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](#), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited



Protocol status: Working

We use this protocol and it's working

Created: December 13, 2022

Last Modified: January 09, 2023

Protocol Integer ID: 73935

Keywords: photosynthetic sea slugs protocol for dnazol, photosynthetic sea slugs protocol, extraction of high molecular weight dna, photosynthetic sea slug, wisecaver lab dnazol, dnazol, high molecular weight dna, high quality of dna, weight dna, dna, oxford nanopore, nanopore, extraction

Abstract


Protocol for DNAzol-based extraction of high molecular weight DNA from photosynthetic sea slugs. This streamlined protocol takes less time compared to the lab's prior CTAB-based protocol while still yielding high quality of DNA for Illumina and Oxford Nanopore sequencing.

Troubleshooting




Prepare Station and Materials

45m

- 1 Set the ThermoMixer to  35 °C
- 2 Get liquid nitrogen.
- 3 Wipe down your work area with 70% ethanol:
 - Chemical fume hood
 - Benchtop
 - Rotors on the microcentrifuge and bench top spinner
 - Pipettes

Prepare slug sample(s)






35m

- 4 Freeze slug tissue in liquid nitrogen and pulverize using a Tissue Lyser II (Qiagen) or a mortar and pestle
- 5 Immediately add the DNazol (see step 9) **or** store in  -80 °C freezer.


It's important to work quickly here as the liquid nitrogen will have lysed many cells already. Material will be highly susceptible to environmental nucleases if allowed to sit at room temperature.



Homogenization and DNA extraction


45m



- 6 Add  300 µL DNazol. Pipet up and down to mix and transfer to a 1.5 mL centrifuge tube.
- 7 Add  5 µL RNase A. Make sure the cap is closed securely, and mix well by vigorously shaking the tube for  00:00:15 .
- 8 Incubate on the ThermoMixer at  500 rpm, 35°C, 00:20:00 . About half way through the incubation, take the tubes out and again mix well by vigorously shaking the tube for  00:00:15 . Place back in the ThermoMixer to complete the incubation step.



8.1 During incubation, prepare an aliquot of 100% ethanol. Prepare  300 μ L per sample.

8.2 During incubation, prepare an ethanol and DNazol wash by mixing  1 mL DNazol with  750 μ L ethanol.

8.3 During incubation, prepare a 75% ethanol wash using nuclease free water. Prepare  300 μ L per sample.

9 Add  300 μ L chloroform to each sample. Make sure the cap is closed securely, and mix well by vigorously shaking the tube for  00:00:15 .


15s




All liquid chloroform liquid waste should be ejected or decanted into the amber bottle labeled 'Phenol:Chloroform Waste' inside the fume hood. Any plasticware that comes into contact with chloroform needs to be properly disposed of in the sharps bin inside the fume hood labeled 'Mutagen/Carcinogen Sharps Do Not Autoclave'.

10 Incubate on the Hula Mixer for  00:05:00

5m

11 Centrifuge at  12000 rpm, 00:10:00 . Don't jostle samples when removing from the centrifuge.


10m

12 Transfer approximately  300 μ L of the top aqueous phase to a fresh 1.5 mL tube. Pipet slowly from the top, being careful not to disrupt the interphase.




DNA Precipitation

35m

13 Add  225 μ L of 100% ethanol. Invert ten times to mix.

14 Incubate at room temperature for  00:05:00 .

5m

15 Centrifuge at  5000 rpm, 00:04:00 . Orient the centrifuge tube so that cap hinge is on the outer edge of the rotor.

4m

Proper orientation of the tube helps to avoid the pelleted RNA when removing the supernatant in subsequent steps. Even if you can't see the pellet, you will know that it is




located on the on the bottom corner under the hinge.

- 16 Discard the supernatant by taking off the initial volume with a 1 mL pipette tip. Use a 200 mL pipette tip to get the last bit of liquid.




This is the last supernatant and plastic waste that needs to be disposed of in the phenol:chloroform waste. All waste in the remaining steps can go in the standard lab 'look-alike' waste bin.

- 17 Add  300 μL of previously prepared DNazol wash and drag along the top of a tube rack to agitate the sample and dislodge the pellet.

- 18 Incubate at room temperature for  00:05:00 .


5m


- 19 Centrifuge at  5000 rpm, 00:04:00 . Orient the centrifuge tube so that cap hinge is on the outer edge of the rotor.

4m

- 20 Discard the supernatant by taking off the initial volume with a 1 mL pipette tip. Use a 200 mL pipette tip to get the last bit of liquid.

Be sure to avoid the hinge side of the tube where the sample is.


- 21 Add  300 μL 75% ethanol wash and vortex briefly. If the pellet does not dislodge at this step, that is okay.

- 22 Centrifuge at  5000 rpm, 00:04:00 . Orient the centrifuge tube so that cap hinge is on the outer edge of the rotor.

4m

- 23 Discard the supernatant by taking off the initial volume with a 1 mL pipette tip. Briefly spin down the tube using the little bench top spinner, and then use a 200 mL pipet tip to get the last bit of liquid.



Again, even if you can't see the pellet, you will know that it is located on the on the bottom corner under the hinge. So be careful to avoid that side of the tube. Pipette away the supernatant by tipping the tube slightly towards you and pipetting the liquid from the bottom corner opposite the hinge.

- 24 Air dry the DNA pellet for  00:05:00 by placing the tube on its side inside a folded kim wipe.

5m



Do not over dry the DNA pellet as this will greatly decrease its solubility. However, also be sure that all the ethanol is removed as residual ethanol will also decrease solubility.

- 25 Resuspend sample in  50 μ L TE Buffer.
- 26 Incubate on the ThermoMixer  500 rpm, 35°C, 00:05:00 .
- 27 Briefly spin down the tube using the little bench top spinner.

QC and clean up

35m

- 28 Make sure all chloroform waste is properly disposed of in the chemical fume hood. Wipe down the chemical fume hood and the benchtop with 70-75% ethanol.
- 29 Assess DNA integrity using the TapeStation. Make sure you reserve time on the TapeStation Teams calendar.
- 30 While the TapeStation is running, measure the DNA concentration using the Qubit.
- 31 If the TapeStation reports a high DNA integrity (DIN) score (≥ 7) and DNA concentration ≥ 20 ng/ μ L, then you have successfully isolated DNA for whole genome sequencing!

Cap your sample tube tightly. Make sure each tube is labeled on the cap and vertically on the side with the sample's unique identifier*. Wrap the cap with parafilm and place sample in the 'Novogene' box in the full sized -20 freezer.

**The identifier should start with a letter and be no more than 12 characters long. Write it carefully using a fine tip permanent marker as this tube will be shipped off for DNA sequencing.*