

Jan 28, 2022

X-ray Micro Computed Tomography (MicroCT)

DOI

<https://dx.doi.org/10.17504/protocols.io.b35aqq2e>



Marta Wawrzyniak¹, Nathalie Weber¹, Simon Blanchoud¹

¹University of Fribourg

Blanchoud lab, UNIFR



Marta Wawrzyniak

University of Fribourg

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.b35aqq2e>

Protocol Citation: Marta Wawrzyniak, Nathalie Weber, Simon Blanchoud 2022. X-ray Micro Computed Tomography (MicroCT) . protocols.io <https://dx.doi.org/10.17504/protocols.io.b35aqq2e>



License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: January 21, 2022

Last Modified: January 28, 2022

Protocol Integer ID: 57218

Keywords: Colonial tunicates, MicroCT, tomography, ascidians, microct for comparative morphology, colonial tunicate, microct, 3d imaging, tissue, simple staining method, micro

Abstract

This protocol has been successfully used with colonial tunicates. It has been adapted to our needs based on the following publication: " MicroCT for comparative morphology: simple staining methods allow high-contrast 3D imaging of diverse non-mineralized animal tissues " by Brian D Metscher.

MicroCT for comparative morphology: simple staining methods allow high-contrast 3D imaging of diverse non-mineralized animal tissues.



Materials

4% PFA

PTA

IKI

Ethanol

low melting agarose

fixing paste (e.g. dental wax)

pipette tips 200uL or 1mL

10x PBS, pH 7.2 (store for up to 2 months at RT)

	A	B	C
	Na ₂ HPO ₄ *2 H ₂ O	0.1M	17.8g
	KH ₂ PO ₄	18mM	2.4g
	NaCl	1.4M	80g
	KCl	27mM	2g
	water		1000mL

10x PBS (amounts calculated for 1L)



Fixation

1h 30m

- 1 Clean the slide with the colony of your interest. See [Cleaning colonial ascidians V.2](#).
- 2 Anesthetize the animals following this protocol. See [Whole colony fixation V.1](#)
- 3 Fix the animals in 4% PFA  Overnight at  Room temperature . See [Whole colony fixation V.1](#)
- 4 Wash twice in 3.3x PBS for  00:30:00 . 30m
- 5 Rinse three times with 1x PBS and leave in 1x PBS for  01:00:00 . 1h

Staining

4d

- 6 Detach the colony with a razor blade and put in a 2mL tube containing  1 mL PTA /  0.6 mL IKI/  0.4 mL Ethanol.
- 7 Incubate for  96:00:00 at  Room temperature on linear shaker (low speed), protect from light. 4d
- 8 Rinse quickly 3-4 times with 1x PBS.

Mounting

- 9 Prepare  10 mL of  0.5 % volume low melting agarose.



- 9.1 When the agarose has cooled down a bit, pipette  125-150 μL if using a 200ul tip or  750-800 μL if using a 1mL tip (the tip size depends on the animal size).
- 9.2 Carefully place your colony vertically inside the tip (it should be as flat as possible, avoid folds).
- 9.3 Pipette some more agarose in order to cover the animal entirely.
- 10 Place the tip into a beaker filled with ice-cubes so that the agarose becomes solid.
- 11 Cover the tip with fixing paste.
- 12 Store at  4 °C or proceed to the next step.
- 13 Mount the tip on the tomograph support using fixing paste in such way that it is perfectly straight.
- 14 Place it inside the tomograph.
- 15 Follow the tomograph's manual and proceed to scan.
- 16 After scanning, the tip with the sample can be stored at  4 °C for several weeks and eventually can be reused.