

Jul 10, 2024

Yeast protoplast fusion

DOI

<https://dx.doi.org/10.17504/protocols.io.5qpvokrxdl4o/v1>

is Sparrow¹, Ulschan Bathe², Kristen Van Gelder²

¹UIUC; ²UF

Yeast Protocols, Tools, a...



is Sparrow

UIUC

Create & collaborate more with a free account

Edit and publish protocols, collaborate in communities, share insights through comments, and track progress with run records.

Create free account

OPEN  ACCESS



DOI: <https://dx.doi.org/10.17504/protocols.io.5qpvokrxdl4o/v1>

Protocol Citation: is Sparrow, Ulschan Bathe, Kristen Van Gelder 2024. Yeast protoplast fusion. **protocols.io**
<https://dx.doi.org/10.17504/protocols.io.5qpvokrxdl4o/v1>

License: This is an open access protocol distributed under the terms of the **[Creative Commons Attribution License](#)**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: In development

We are still developing and optimizing this protocol

Created: May 13, 2024

Last Modified: July 10, 2024

Protocol Integer ID: 99664

Keywords: yeast protoplast fusion this protocol, yeast protoplast fusion, yeast protoplast, yeast strain, yeast, fusion, hanson lab, original protocol by ulschan bathe, original protocol

Disclaimer

DISCLAIMER – FOR INFORMATIONAL PURPOSES ONLY; USE AT YOUR OWN RISK

The protocol content here is for informational purposes only and does not constitute legal, medical, clinical, or safety advice, or otherwise; content added to protocols.io is not peer reviewed and may not have undergone a formal approval of any kind. Information presented in this protocol should not substitute for independent professional judgment, advice, diagnosis, or treatment. Any action you take or refrain from taking using or relying upon the information presented here is strictly at your own risk. You agree that neither the Company nor any of the authors, contributors, administrators, or anyone else associated with protocols.io, can be held responsible for your use of the information contained in or linked to this protocol or any of our Sites/Apps and Services.

Abstract

This protocol was based on the original protocol by Ulschan Bathe and Kristen Van Gelder from the Hanson Lab at UF in 2023.

It details how to fuse 2 yeast protoplasts. It is focused on obtaining an evolution strain with the system OrthoRep, but should work for the fusion of any 2 yeast strains provided the selection scheme is functional.



Media and buffers

- 1 Make the following solutions ahead of time. They do not need to be sterilized.

Unless specified, all solutions are made in water

1.1 CPB (Citrate Phosphate Buffer) solution

Solution A

[M] 0.1 Mass Percent Citric Acid (dihydrate)

Solution B

[M] 0.2 Mass Percent Na_2HPO_4

Combine  34.8 mL A and  65.2 mL B and bring up final volume to  200 mL with MilliQ water

1.2 3M KCl

1.3 0.5 M EDTA (pH = 8.0)

1.4 0.5 M CaCl_2

1.5 Other media

YPD

Autoclaved sterile water

- 2 Solutions - made fresh the day of

2.1 BMEE solution

 3 mL [M] 0.5 Molarity (M) EDTA

 50 μL beta-mercaptoethanol



 21.95 mL MilliQ water

Filter sterilize

2.2 0.6 M KCl

 20 mL  3 Mass Percent KCl

 80 mL MilliQ water

Filter sterilize

2.3 Buffer 1

 39 mL CPB Buffer

 10 mL  3 Mass Percent KCl

 1 mL  0.5 Mass Percent EDTA

Filter sterilize

2.4 Buffer 2

A -  50 % w/v **PEG 3350**

Add  10 g PEG +  8 mL water, stir with high heat and adjust volume to 25 mL when done

B - Buffer 2

 16.65 mL  50 % w/v PEG 3350

 5 mL  3 Mass Percent KCl

 2.5 mL  0.5 Mass Percent CaCl₂

 850 µL MilliQ water

Filter sterilize

2.5 Buffer 3

 5 mL  3 Mass Percent KCl

 2.5 mL  0.5 Mass Percent CaCl₂



 17.5 mL MilliQ water

Filter sterilize

Plating media

- 3 Plating media (i.e., selective media) must be made in 100-150 mL batches. Each 25 mL supports a single plating, thus 100 mL supports 1 fusion with 2 platings at different concentrations

For  100 mL

 93.5 mL MilliQ water

 4.473 g **KCl**

 0.667 g YNB with ammonium sulfate OR  0.5 g ammonium sulfate +  0.171 g

YNB without ammonium sulfate (monosodium glutamate can also be used as an alternative nitrogen source if ammonium sulfate cannot be used due to interfering with antibiotics (G418 or Nat))

 0.14 g g of appropriate dropout

Adjust pH to 6.1

 3 g agar

Run on liquid 30 cycle, leave inside autoclave until later.

Yeast culture

- 4 2 or 3 days before the experiment, start  3 mL pre-cultures in appropriate selection media for the donor (**p1 carrying strain**) and recipient (**background + EP-DNAP**) strains. Grow at  30 °C with shaking

2d

- 5 The day before protoplast fusion, dilute pre-cultures 1:50 into  50 mL YPD culture and incubate shaking at  30 °C

16h

Be very mindful of the rate of growth of the cells - OD after outgrowth should not exceed 7 but should be above 2. Depending on your strain, you will need to figure out optimal time for starting the culture.

Protoplast fusion

1w 3d 3h 33m

- 6 The morning of protoplast fusion, make the solutions above and filter-sterilize them.



- 7 Take  100 μ L of culture and add  900 μ L YPD to it, then measure the OD600.

Multiply the OD by the dilution factor (x10)

Using the approximation that, at OD600 = 4.5, for 20 mL of culture, ~0.45 g of cells are yielded, calculate the volume of culture necessary to obtain  0.3 g of cells of each strain using the formula **$c1.v1 = c2.v2$**

- 8 Weigh your tubes before adding culture and make a note of it.

- 9 Centrifuge at  3000 x g, 00:05:00 , using blank tubes to balance the different culture volumes

5m

- 10 Weigh your tubes again, and adjust pellet weight by adding more culture or scooping out pellet.

Aim for a final weight of  0.3 g - preferably OVER

- 11 Resuspend in 10 mL sterile water and transfer to 15 mL falcon

- 12 Centrifuge at  3000 x g, 00:05:00 , and decant supernatant. Pipette out water.

5m

- 13 Resuspend in  1.8 mL in the same tube **BMEE solution TAPPING GENTLY to resuspend**

DO NOT VORTEX ANYTHING BEYOND THIS POINT

- 14 Incubate at  30 $^{\circ}$ C for  00:30:00 .

30m

At 15 min, invert a few times to mix gently.

- 15 Centrifuge at  3000 x g, 00:05:00 , decant supernatant.

5m



16 Add  3 mL **0.6 M KCl** and resuspend with serological pipette.

Note: avoid the creation of bubbles

17 Centrifuge at  3000 x g, 00:05:00 , decant supernatant.

5m

18 Add  4.8 mL **Buffer I** and resuspend with serological pipette.

19 Add Zymolyase solution at a total concentration of  6 U/mL

Note: Hanson Lab storage buffer for Zymolyase is as follows. If using this recipe make at a concentration of 5 U/uL, then add 5.8 uL :

- i. 100 mg (2000 U) of zymoylase from amsbio (Cat# = 120491-1)
- ii. 0.1 M Na₂PO₄ (pH 7.5, adjusted with NaOH)
- iii. 50% glycerol
- iv. 3 mM β-ME

To make 400 uL of solution

1. 100 mg zymolyase
2. 200 uL glycerol
3. 192 uL 0.1 M Na₂PO₄ (pH 7.5, adjusted with NaOH)
4. 8 uL of a 1:100 beta-mercaptoethanol dilution in 0.1 M Na₂PO₄

20 Incubate at  30 °C statically for  01:00:00 and rotate every  00:15:00 by hand.

1h 15m

Use these incubations to make the plating media and autoclave it.

Additionally, set up your  42 °C water bath next to the fume hood. Use a thermometer to make sure the temperature is right.

21 Centrifuge at  700 x g, 0°C, 00:10:00 , decant supernatant.

30m



Add  3 mL **0.6 M KCl**. Resuspend gently by adding the buffer along the side of the tube.

Centrifuge at  700 x g, 0°C, 00:10:00 , decant supernatant.

Add  3 mL **0.6 M KCl**. Resuspend gently by adding the buffer along the side of the tube.

Centrifuge at  700 x g, 0°C, 00:10:00 , decant supernatant.

22 Resuspend in  3 mL **Buffer 1**

23 In new 15 mL tubes, mix  1.5 mL of donor strain with  1.5 mL of recipient strain.

	A	B	C
		Donor strain	Donor strain
	Recipient strain	Evolution strain	Evolution strain
	NA	Control 1	-

Use appropriate EP-DNAPs

The remaining volumes can be carried forward as negative controls.

Mix by inversion gently a few times.

24 Centrifuge at  700 x g, 0°C, 00:10:00 , decant supernatant. Resuspend in  5 mL of Buffer 2.

10m

The PEG will make resuspension difficult, so try your best with the serological pipette. Tiny, visible clumps of cells are OK, but should be avoided as they can make identification of emerging colonies difficult later on - avoid using the small pipettes.

25 Incubate at  30 °C for  00:30:00 with occasional inverting.

30m



26 Centrifuge at  700 x g, 0°C, 00:10:00 , decant supernatant.

15m

The media should be done autoclaving at this point. Take out, add glucose and other required auxotrophic additives, and return back to the autoclave to maintain liquidity for at least  00:05:00

BE VERY CAREFUL NOT TO ADD ADDITIVES WHICH LOSE YOUR SELECTION SCHEME!!!! - PAY ATTENTION TO SELECTION MARKER IN P1 (LEU2) AND SELECTION MARKER IN EP-DNAP CONSTRUCT (In our case, this is URA3, but other plasmids have HIS3)

27 Resuspend in  5 mL **Buffer 3** with EXTREME GENTLENESS - the cells are protoplasted at this point.

28 Take media in the autoclave and place in  42 °C water bath, for a minimum of  00:03:00 . **TIME IT!!!!!!**

3m

You should be able to touch the flask with your wrist's bare skin and not feel pain, only warmth.

29 In a 50 mL falcon tube, add cells

1w 3d

Make 2-3 tubes per fusion:

- One with  250 µL of cells
- One with  500 µL of cells
- One with  1 mL of cells

Gently, add 25 mL of plating media, mix gently, and then pour in labelled plates. Allow plates to solidify and incubate at  30 °C for up to  240:00:00

30 On the day after, tape the lid with parafilm.