



Aug 16, 2019

Zona pellucida removal using Acid Tyrode solution and snap freezing

DOI

dx.doi.org/10.17504/protocols.io.6kwhcxe



Elizabeth Jannaman¹, Peter J Hansen¹

¹University of Florida

University of Florida Bovine Embryo Lab
Tech. support email: hansen@animal.ufl.edu



Peter J Hansen

University of Florida

OPEN  ACCESS



DOI: dx.doi.org/10.17504/protocols.io.6kwhcxe

External link: https://animal.ifas.ufl.edu/hansen/lab_protocol_docs/Acidtyrodes.pdf

Protocol Citation: Elizabeth Jannaman, Peter J Hansen 2019. Zona pellucida removal using Acid Tyrode solution and snap freezing. **protocols.io** <https://dx.doi.org/10.17504/protocols.io.6kwhcxe>

License: This is an open access protocol distributed under the terms of the **Creative Commons Attribution License**, which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited

Protocol status: Working

We use this protocol and it's working

Created: August 16, 2019

Last Modified: August 16, 2019

Protocol Integer ID: 26998

Keywords: preimplantation embryo, bovine, zona pellucida, Acid Tyrode's



Abstract

A simple protocol to remove the zona pellucida from bovine embryos. This is a necessary step before freezing embryos for analysis of RNA or DNA because of contaminating cumulus and sperm cells associated with the zona pellucida.

Attachments



[Acidtyrodes.pdf](#)

90KB

1